

M5 HiPure pET21a(+)表达载体/vector, C-末端 His6 标签 使用说明书

Product	Unit	Cat.#
M5 HiPure pET21a(+)表达载体/vector	10 µl, 20ng/ul	MFV01001-01


保存: -20℃ 保存

描述: C-末端 His6 标签

成分: 10 mM Tris (pH8.0)

序列信息:

pET21a(+)

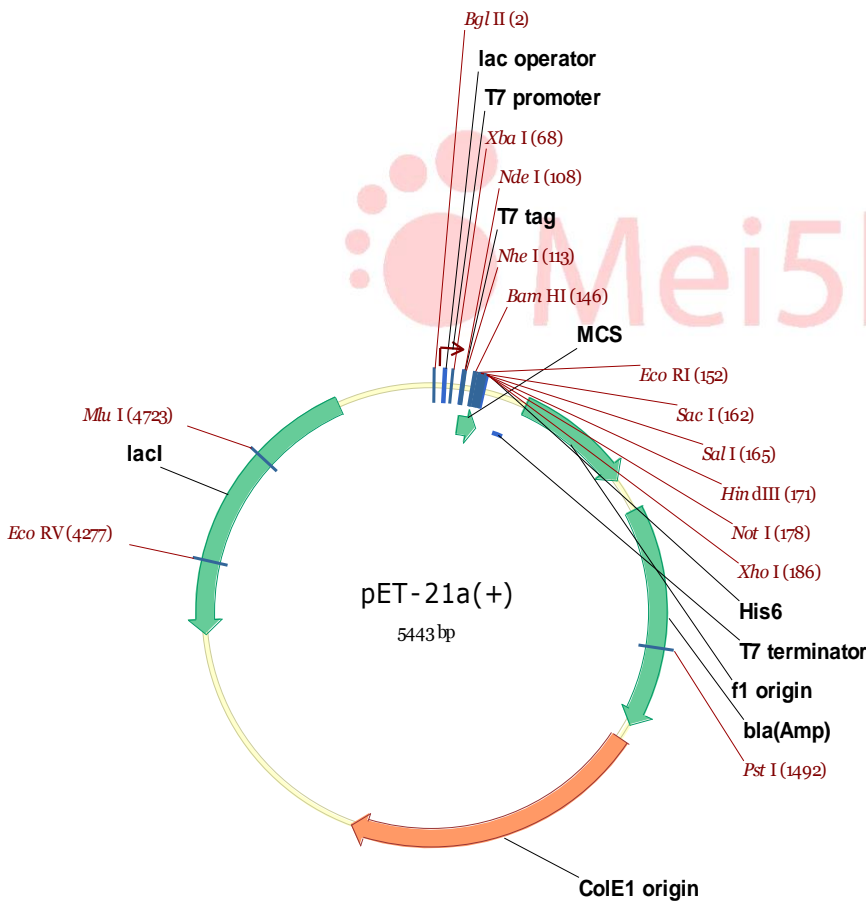

LOCUS pET-21a(+) 5443 bp DNA circular 23-JUL-2013
 SOURCE
 ORGANISM
 COMMENT This file is created by Vector NTI
 COMMENT ORIGDB|GenBank

```

COMMENT      VNTDBDATE|668342873|
COMMENT      LSOWNER|
COMMENT      VNTNAME|pET-21a(+)|
FEATURES      Location/Qualifiers
    misc_feature   946..1803
                  /vntifkey="21"
                  /label=bla(Amp)
    misc_feature   359..814
                  /vntifkey="21"
                  /label=f1\origin
    promoter       21..37
                  /vntifkey="30"
                  /label=T7\promoter
    misc_feature   40..64
                  /vntifkey="21"
                  /label=lac\operator
    terminator     276..322
                  /vntifkey="43"
                  /label=T7\terminator
    misc_feature   106..190
                  /vntifkey="21"
                  /label=MCS
    misc_feature   191..208
                  /vntifkey="21"
                  /label=His6
    misc_feature   complement(3998..5077)
                  /vntifkey="21"
                  /label=lacI
    misc_feature   109..141
                  /vntifkey="21"
                  /label=T7>tag
    rep_origin     1867..3034
                  /vntifkey="33"
                  /label=ColE1\origin
BASE COUNT    1268 a      1441 c      1446 g      1288 t
ORIGIN

```

```

1 agatctcgat cccgcgaaat taatacact cactataggg gaattgtgag cggataacaa
61 tcccctcta gaaataatt tgttaactt taagaaggag atatacatat ggctagcatg
121 actggtggac agcaaatggg tcgcggatcc gaattcgagc tccgtcgaca agcttgcggc
181 cgcactcgag caccaccacc accaccactg agatccggct gctaacaaag cccgaaagga
241 agctgagttg gctgctgcca ccgctgagca ataactagca taacccttg gggcctctaa
301 acgggtcttg aggggtttt tgctgaaagg aggaactata tccggattgg cgaatgggac
361 gcgccctgta gggcgcatt aagcgcggcg ggtgtggtgg ttacgcgcag cgtgaccgct
421 acactgcca ggccttagc gcccgctct ttcgcttct tccttctt tctcgccacg
481 ttcgccggt tcccctgca agctctaaat cgggggctcc ctttaggggt cegatttagt
541 gctttacggc acctcgacc caaaaaactt gattagggtg atggttcacg tagtgggcca
601 tcgccctgat agacggttt tcgcccttg acgttggagt ccacgttctt taatagtgga
661 ctctgttcc aaactggaac aacctcaac cctatctcgg tctattctt tgattataa
721 gggattttgc cgatttcggc ctattggtta aaaatgagc tgatttaaca aaaattaac
781 gcgaatttta acaaaatatt aacgtttaca attcaggtg gcacttttcg gggaaatgtg
841 cgcggaacce ctattgttt atttttctaa atacattcaa atatgtatcc gctcatgaga
901 caataacctt gataaatgct tcaataatat tgaaaaagga agagtatgag tattcaacat
961 ttcggtgag cccttatcc ctttttgcg gcattttgcc ttctgtttt tgcaccaca
1021 gaaacgctgg tgaagtaaaa agatgctgaa gatcagttgg gtgcacgagt gggttacatc
1081 gaactggatc tcaacagcgg taagatcctt gagagtttc gccccgaaga acgtttcca

```

1141 atgatgagca cttttaaagt tctgctatgt ggcgcggtat tatcccgtat tgacgccggg
1201 caagagcaac tcggtcgcgc catacactat tctcagaatg acttggttga gtactacca
1261 gtcacagaaa agcatcttac ggatggcatg acagtaagag aattatgcag tgctgccata
1321 accatgagtg ataactctgc ggccaactta ctctgacaa cgatcggagg accgaaggag
1381 ctaaccgctt tttgcacaa catgggggat catgtaactc gccttgatcg ttgggaaccg
1441 gagctgaatg aagccatacc aaacgacgag cgtgacacca cgatgcctgc agcaatggca
1501 acaacgttgc gcaaactatt aactggcga a ctacttact tagctccc gcaacaatta
1561 atagactgga tggaggcggg taaagttgca ggaccactc tgcgctcggc ccttcggct
1621 ggctggttta ttgctgataa atctggagcc ggtgagcgtg ggtctcgcgg tatcattgca
1681 gcaactatgg atgaacgaaa tagacagatc gctgagatag gtcctcact gattaagcat
1741 gcaactatgg atgaacgaaa tagacagatc gctgagatag gtcctcact gattaagcat
1801 tggtaactgt cagaccaagt ttactcatat atactttaga ttgattaaa acttcattt
1861 taatttaaaa gtagctaggt gaagatcctt ttgataatc tcatgacca aatcccttaa
1921 cgtgagtttt cgttccactg agcgtcagac cccgtagaaa agatcaaagg atcttctga
1981 gatccttttt ttctgcgctg aatctgctgc ttgcaaacaa aaaaaccacc gctaccagc
2041 gtggtttgtt tgcccggata agagctacca actcttttc cgaaggtaac tggcttcagc
2101 agagcgcaga taccaaatc tgccttcta gtgtagccgt agttaggcca ccaactcaag
2161 aactctgtag caccgcctac atacctcgtc ctgtaatcc tgttaccagt ggctgctgcc
2221 agtggcgata agtcgtgtct taccgggtg gactcaagac gatagttacc ggataaggcg
2281 cagcggtcgg gctgaacggg gggtctgctc acacagccca gcttggagcg aacacctac
2341 accgaactga gatactaca gcgtgagcta tgagaaagcg ccacgctcc cgaagggaga
2401 aagcgggaca ggtatccggt aagcggcagg gtcggaacag gagagcgcac gagggagctt
2461 ccagggggaa acgcttgta tctttatagt cctgctgggt ttcgccacct ctgactgag
2521 cgtcgtatft tctgatgctc gtcagggggg cggagcctat ggaaaaacgc cagcaacgc
2581 gccttttac ggttctgctc cttttgctg cttttgctc acatgtctt tctgctgta
2641 tcccctgatt ctgtggataa ccgtattacc gcctttgagt gagctgatac cgtcgcgc
2701 agccgaacga ccgagcgcag cgagtcagt agcaggaag cgaagagcg cctgatgcgg
2761 tattttctcc ttacgcatc gtgcggtatt tcacaccga tatatggtgc acttcagta
2821 caatctgctc tgatccgca tagttaagcc agtatacact ccgctatgc tacgtgactg
2881 ggtcatggtt gcgccccgac acccgccaac acccgtgac gcgccctgac gggcttctc
2941 gctcccggca tccgcttaca gacaagctgt gaccgtctc gggagctgca tctgtcagag
3001 gttttaccg tcataccga aacgcgcgag gcagctcgg taaagctcat cagcgtgctc
3061 gtgaagcgat tcacagatgt ctgcctgttc atccgcgtcc agctcgttga gttctccag
3121 aagcgttaat gctcggctc tgataaagcg ggccatgta agggcgggtt ttctctgtt
3181 ggtcactgat gcctcctgtt aagggggatt tctgttcag ggggtaatga taccgatgaa
3241 acgagagagg atgctcacga tacgggttac tgatgatgaa catgcccggt tactggaacg
3301 ttgtgagggt aaacaactgg cggtatggt ggcggcggac cagagaaaa tcaactcagg
3361 tcaatgccag cgtctgta atacagatgt aggtgtcca cagggtagcc agcagcatcc
3421 tgcgatgcag atccggaaca taatggtgca gggcgtgac ttccgcttt ccagactta
3481 cgaaacacgg aaaccgaaga ccattcatgt tttgtctag gtcgcagacg tttgcagca
3541 gcagtcgctt cagcttctc cgcgtatcgg tgattcattc tgctaaccag taaggcaacc
3601 ccgccagcct agccgggtcc tcaacgacag gagcacgac atgcgcacc gtggggccgc
3661 catccggcg ataatggcct gttctcgc gaaacgtttg gtggcgggac cagtgcgaa
3721 ggcttgagcg agggcgtgca agattccgaa taccgcaagc gacaggccga tcatctcgc
3781 gctccagcga aagcgtctc cccgaaaat gaccagagc gctgccgca cctgtctac
3841 gattgcatg ataaagaaga cagtcataag tgcggcgacg atagtcatgc cccgcgcca
3901 ccggaaggag ctgactgggt tgaaggctct caaggcctc ggtcagatc ccggtgcta
3961 atgagtgagc taactacat taattgctt gcgctcact cccgcttcc agtcgggaaa
4021 cctgtcgtc cagctcatt aatgaatcgg ccaacgcgcg gggagaggcg gtttgcgat
4081 tggcgccag ggtggtttt cttttacca gtgagacggg caacagctga ttgccctca
4141 ccgctggcc ctgagagatg tgcagcaagc ggtccacgct ggtttgccc agcaggcgaa
4201 aatcctgttt gatgtggtt aacggcggga tataacatga gctgtcttc gtatcgtct
4261 atcccactc cgagatattc gcaccaacgc gcagcccgga ctcggtaatg gcgcgcttg
4321 cggccagcgc catctgatc ttggcaacca gcatcgcagt gggaacgatg ccctcattc
4381 gcatttgcatt gttttgttga aaaccggaca tggcactcca gtcgcttcc cgttccgcta

4441 tcggctgaat ttgattgca gtgagatatt tatgccagcc agccagacgc agacgcgccg
4501 agacagaact taatgggcc gctaacacgc cgatttctg gtgaccaat gcgaccagat
4561 gctccagcc cagtcgcgta ccgttctcat gggagaaaat aatactgtg atgggtgtct
4621 ggtcagagac atcaagaat aacgccgga cattagtca ggcagctcc acagcaatgg
4681 catctggtc atccagcga tagttaatga tcagcccact gacgcgtgc gcgagaagat
4741 tgtcaccgc cgcttacag gcttcgacgc cgcttcgtc taccatgac accaccacgc
4801 tggcaccag ttgatcgccg cgagattaa tcgccgcgac aatttgcgac ggcgcgtgca
4861 gggccagact ggaggtggca acgcaatca gcaacgactg tttgccgcc agttgtgtg
4921 ccacgcggtt gggaatgta ttcagctccg ccacgcgcgc tccacttt tcccgcgtt
4981 tcgcaaaac gtggctggcc tggttacca cgcgggaaac ggtctgataa gagacaccgg
5041 catactctg gacatcgtat aacgttactg gttcacatt caccacctg aattgactct
5101 cttccggcg ctatcatgcc ataccgcga aggttttgcg ccattcgatg gtgtccggga
5161 tctcgacgt ctccttatg cgactcctgc attaggaagc agcccagtag taggtgagg
5221 cggttgagca ccgccgccg aaggaatgt gcatgcaagg agatgggcc caacagtccc
5281 ccggccacgg ggctgccac catacccac cgaaacaag cgctcatgag cccgaagtgg
5341 cgagccgat ctcccctc ggtgatgctg gcgatatagg cgccagcaac cgcacctgtg
5401 gcgccgtga tgccggccac gatgcgtccg gcgtagagga tcg

//



Please note: All products are "FOR RESEARCH USE ONLY AND ARE NOT INTENDED FOR DIAGNOSTIC OR THERAPEUTIC USE"

