

---

## Product Information

### **CONCANAVALIN A** **Mei5bio Prod. No. MF104**

**CAS NUMBER:** 11028-71-0

**SYNONYM:** Con-A

#### **PHYSICAL PROPERTIES:**

Appearance: white to white with a yellow cast powder

Molecular Weight: 25,500<sup>1</sup> (see Description section for additional details)

Extinction Coefficient:  $E^{1\%}(280\text{nm}) = 11.4$  (0.1 M NaCl)<sup>2</sup>

$E^{1\%}(280\text{nm}) = 13.7$  (0.05 M sodium phosphate pH 6.8 with 0.2 M NaCl)

$E^{1\%}(280\text{nm}) = 12.4$  (0.05 M sodium acetate pH 5.2 with 0.2 M NaCl)<sup>3</sup>

Isoelectric point:  $pI = 4.5, 4.7, 5.0-5.1, \text{ and } 5.4-5.5$  (several isoforms)<sup>4</sup>

#### **DESCRIPTION:**

Con-A is a lectin isolated from the jack bean. Con-A is not a glycoprotein.<sup>5</sup> The monomeric molecular weight of Con-A is 25,500.<sup>1</sup> At pH 5.5 Con-A exists as a dimer, and at pH>7 it exists as a tetramer.<sup>6</sup> Con-A does not contain cysteine residues.<sup>2</sup> Unlike most other lectins, Con-A is a metalloprotein and requires a transition metal ion, such as manganese, plus calcium ions for binding.<sup>7</sup>

#### **STABILITY / STORAGE AS SUPPLIED:**

Con-A has a shelf-life of three years when stored properly below 0°C. At Mei5bio agglutination titer of a three-year-old lot was identical to the original assay value.

#### **SOLUBILITY / SOLUTION STABILITY:**

Mei5bio tests the solubility of MF104 at 10 mg/ml in water which produces a slightly hazy to hazy colorless to faint yellow solution. The solubility of Con-A improves with increasing concentrations of sodium chloride up to 1 molar.

## CONCANAVALIN A Mei5bio Prod. No. MF104

### SOLUBILITY / SOLUTION STABILITY: (continued)

Solution stability is also improved with increasing concentrations of sodium chloride. The following solution stability data was obtained using a 20 mM Tris HCl buffer with 1 mM CaCl<sub>2</sub> plus 0.5 mM MnCl<sub>2</sub> pH 7.0 with the stated concentrations of sodium chloride. Less than 5% loss of activity was observed after 10 days at room temperature in buffer containing 1 M NaCl. An 8% loss in activity was observed after 10 days at room temperature in buffer containing 0.2 M NaCl. A 15% loss in activity was observed after 10 days at room temperature in buffer containing 0.1 M NaCl. Zero percent loss in activity was observed after four to six weeks at 0-4 °C in a 0.1 M Tris + 80 mM Glycine buffer pH 5.0 with 3 mM CaCl<sub>2</sub> and 3 mM MnCl<sub>2</sub>.<sup>8</sup>

Solutions of Con A in 1 M NaCl can be frozen and thawed without appreciable loss of activity. KCl should NOT be substituted for NaCl.<sup>10</sup>

Con A has been reported to withstand heat treatment at 45 °C for two hours at pH 3.0-3.2 and at 6-8 °C for 16-18 hours.<sup>9</sup>

### APPLICATIONS:

Con-A binds specifically to mannosyl and glucosyl residues of polysaccharides and glycoproteins.<sup>11</sup> Unmodified hydroxyl groups at the C3, C4 and C6 positions of D-glucopyranosyl or D-mannopyranosyl rings may be essential for binding.<sup>12</sup> Each subunit of Con-A contains one calcium ion and one manganese ion. Removal of these cations by dialysis under acidic conditions abolishes the carbohydrate-binding activity.<sup>11</sup>

Con-A has been used to elucidate structural changes in the membrane surface of transformed cells.<sup>13,14,15,16</sup>

Con-A has been found to have mitogenic activity.<sup>18</sup>

Con-A conjugates of antitumor drugs have been used in drug delivery systems in cultured cells.<sup>19,20,21</sup>

### REFERENCES

1. Wang, J. L., et al., *J. Biol. Chem.*, 250, 1503 (1975).
2. Agrawal, B. and Goldstein, I.J., *Arch. Biochem. Biophys.*, 124, 218 (1968).
3. Yariv, j., et al., *Biochim. Biophys. Acta*, 165, 303 (1968).
4. Entlicher, J.V., et al., *Biochim. Biophys. Acta*, 236, 795 (1971).
5. Olson, M.O. and Liener, I.E., *Biochemistry*, 6, 105 (1967).
6. Senear, D.F. and Teller, D.C., *Biochemistry*, 20, 3076 (1981).
7. Sumner, J.B. and Howell, S.F., *J. Biol. Chem.*, 115, 583 (1936).
8. Sophianopoulos, A.J. and Sophianopoulos, J.A., *Prep. Biochem.*, 11, 413 (1981).
9. Sigma data
10. *The Lectins*, ed Liener, et al., p. 53 (1986).

**CONCANAVALIN A**  
**Mei5bio Prod. No. MF104**

**REFERENCES:** (continued)

11. Goldstein, I.J., et al., *Biochemistry*, 4, 876 (1965).
12. Yahara, I. and Edelman, G.M., *Proc. Nat. Acad. Sci. USA*, 69, 608 (1972).
13. Shoham, J., et al., *Nature*, 227, 1244 (1970).
14. Inbar, M. and Sachs, L., *Proc. Nat. Acad. Sci. USA*, 63, 1418 (1969).
15. Inbar, M., and Sachs, L., *Nature*, 223, 710 (1969).
16. Burger, M. and Noonan, K.D., *Nature*, 228, 512 (1970).
17. Lin, J., et al., *J. Nat. Cancer Inst.*, 66, 523 (1977).
18. Kitao, T. and Hattori, K., *Nature*, 265, 81 (1977).
19. Tsuruo, T., et al., *Int. J. Cancer*, 26, 81 (1977).

\* In addition to MF104 Mei5bio also offers other grades of Con-A, including tissue culture-tested-sterile Con-A, FITC, TRITC, biotin, peroxidase, and ferritin labeled Con-A, plus a complete line of Con-A-based affinity resins. See the Lectin and Cell Culture sections of the Sigma catalog for complete listings.

**Mei5bio warrants that its products conform to the information contained in this and other Mei5bio publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.**