

M5 HiPure pGEX-4T-1 表达载体 vector, N-末端 GST 融合标签 使用说明书

Product	Unit	Cat.#
M5 HiPure pGEX-4T-1 表达载体 vector	10 µl, 20ng/ul	MFV01022-01

保存: -20°C保存

别名: pGEX-4T-1, pGEX4T1, pGEX 4T 1

质粒类型: 大肠杆菌表达载体

表达水平: 高拷贝

启动子: Tac

克隆方法: 多克隆位点, 限制性内切酶

载体大小: 4969 bp (查看载体序列)

5' 测序引物及序列: pGEX5': GGGCTGGCAAGCCACGTTTGGTG

3' 测序引物及序列: pGEX3': CCGGGAGCTGCATGTGTCAGAGG

载体标签: N-GST

载体抗性: Ampicillin (氨苄青霉素)

备注: 复制子是 pMB1

成分: 10 mM Tris (pH8.0)


Gel electrophoresis of QC control:



1% Agarose Gel QC

Lane 1: pGEX-4T-1; M1: DNA marker DL2000

序列信息:

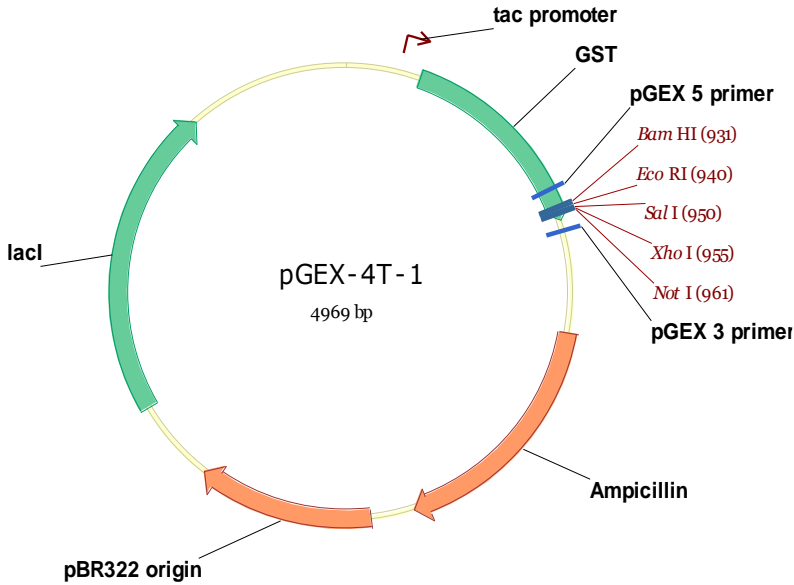

LOCUS	pGEX-4T-1	4969 bp	DNA	circular SYN 16-NOV-2010
DEFINITION	pGEX-4T1.			
SOURCE	ORGANISM other sequences; artificial sequences; vectors.			
FEATURES	Location/Qualifiers			
source	1..4969 /organism="pGEX-4T1" /mol_type="other DNA" /vntifkey="98"			
promoter	184..212 /vntifkey="30" /label=tac_promoter			
misc_feature	258..977 /vntifkey="21" /label=GST			
misc_feature	869..891 /vntifkey="21" /label=pGEX_5_primer			
misc_feature	complement(1019..1041) /vntifkey="21" /label=pGEX_3_primer			
CDS	1377..2237 /gene="Ampicillin" /vntifkey="4" /label=Ampicillin /note="ORF frame 3"			
gene	1377..2237 /gene="Ampicillin" /vntifkey="60"			
rep_origin	2392..3011 /vntifkey="33" /label=pBR322_origin			

```
misc_feature      3309..4400
                  /vntifkey="21"
                  /label=lacI
BASE COUNT      1225 a      1202 c      1292 g      1250 t
ORIGIN
1 acgttatcga ctgcacggcg caccaatgct tctggcgtca ggcagccatc ggaagctgtg
61 gtagtgctgt gcaggtcgtg aatcactgca taatcgtgt cgctcaaggc gcactcccgt
121 tctggataat gtttttgcg cggacatcat aacggttctg gcaaatattc tgaatgagc
181 tgttgacaat taatcatcgg ctctgataat gtgtggaatt gtgagcggat aacaatttca
241 cacaggaaac agtattcatg tcccctatac taggtfattg gaaaattaag ggccttgtgc
301 aacctactcg acttcttttg gaatatcttg aagaaaaata tgaagagcat ttgatgagc
361 gcgatgaagg tgataaatgg cgaacaacaaa agtttgaatt gggtttgag tttccaatc
421 ttcttatta tattgatgtg gatgttaaat taacacagtc tatggccatc atacgttata
481 tagctgaaa gcacaacatg ttgggtggtt gtccaaaaga gcgtgcagag atttcaatgc
541 tgaaggagc ggttttggat attagatacg gtgttcgag aattgcatat agtaaagact
601 tgaactct caaagttgat ttcttagca agctacatg aatgctgaaa atgtcgaag
661 atcgttatg tcataaaaac tatttaaatg gtgatcatgt aacctatcct gacttcatgt
721 tgtatgacg tcttgatgtt gttttataca tggaccaat gtgcctggat gcgttccaa
781 aattagttt ttttaaaaaa cgtattgaag ctatcccaca aattgataag tacttgaat
841 ccagcaagta tatagcatgg ctttgcagg gctggcaagc cacgtttgtt ggtggcgacc
901 atctccaaa atcggatctg gtccgcgctg gatccccgga attccccggt cgactcgagc
961 ggccgcatcg tgaactactg acgatctgcc tcgcgcgttt cggatgatgac ggtgaaaacc
1021 tctgacacat gcagctcccg gagacgggtc cagcttctct gtaagcggat gccgggagca
1081 gacaagcccg tcagggcgcg tcagcgggtg ttggcgggtg tcggggcgca gccatgacc
1141 agtcacgtag cgatagcggg gtgtataat ctgaaagacg aaaggcctc gtgatacgc
1201 tattttata ggttaatgct atgataataa tggtttcta gacgtcaggt ggcactttc
1261 ggggaaatgt gcgcggaacc cctattgtt tattttcta aatacattca aatatgtatc
1321 cgctcatgag acaataacc tgataaatgc ttcaataata tgaaaaagg aagagtatga
1381 gtattcaaca ttccgtgtc gcccttattc cttttttgc ggcattttgc cttctgttt
1441 ttgctcacc agaaacgctg gtgaaagtaa aagatgctga agatcagttg ggtgcacgag
1501 tgggttacat cgaactggat ctcaacgctg gtaagatcct tgagagtttt cgccccgaag
1561 aacgtttcc aatgatgagc acttttaaag ttctgctatg tggcgcggta ttatccgtg
1621 ttgacgcgg gcaagagcaa ctccgtcgc gcatacacta ttctcagaat gacttggtg
1681 agtactcacc agtcacagaa aagcatctta cggatggcat gacagtaaga gaattatgca
1741 gtgctgcat aacctagatg gataacactg cggccaactt actctgaca acgatcggag
1801 gaccgaagga gtaaccgct ttttgcaca acatggggga tcatgtaact cgccttgatc
1861 gttgggaacc ggagctgaa gaagccatac caaacgacga gcgtgacacc acgatgcctg
1921 cagcaatggc aacaacgttg cgcacaactat taactggcga actacttact ctagtccc
1981 ggcaacaatt aatagactgg atggaggcgg ataaagtgc aggaccactt ctgcgctcgg
2041 ccttccggc tggctggtt attgctgata aatctggagc cggtagcgtt gggctcgcg
2101 gtatcattg agcactgggg ccagatggta agccctccc tatcgtatgt atctacacga
2161 cggggagtca ggcaactatg gatgaacgaa atagacagat cgctgagata ggtgcctcac
2221 tgattaagca ttgtaactg tcagaccaag ttactcata tatactttag attgattaa
2281 aactcattt taafttaaa aggatctagg tgaagatcct tttgataat ctatgacca
2341 aaatcccta acgtgagttt tcgtccact gagcgtcaga ccccgtagaa aagatcaaag
2401 gatcttctg agatcctttt ttctgcgcg taatctgctg cttgcaaaac aaaaaccac
2461 cgctaccagc ggtggtttg ttgccgcatc aagagctacc aactctttt ccgaaggtaa
2521 ctggttcag cagagcgcag ataccaata ctgtcctct agttagcgg tagttaggcc
2581 accactcaa gaactctgta gcaccgccta catacctcgc tctgctaact ctgttaccag
2641 tggctgctg cagtggcgat aagtcgtgtc ttaccgggtt ggactcaaga cgatagttac
2701 cggataaggc gcagcggctg ggctgaacgg ggggttcgtg cacacagccc agcttgagc
2761 gaacgaccta caccgaactg agatacctac agcgtgagct atgagaaaac gccacgctc
2821 ccgaaggag aaaggcggac aggtatccgg taagcggcag ggtcggaaac ggagagcgca
2881 cgaggagct tccaggggga aacgcctggt atctttatg tctgtcggg tttcgccacc
2941 tctgactga gcgtcattt ttgtatgct cgtcaggggg gcggagccta tggaaaaacg
```

3001 ccagcaacgc ggccttttta cggttcctgg cttttgctg gcctttgct cacatgtct
3061 ttctgcgtt atcccctgat tctgtgata accgtattac cgcctttgag tgagtgata
3121 ccgctgcgcg cagccgaacg accgagcgca gcgagtcagt gagcgaggaa gcggaagagc
3181 gcctgatgcg gtattttctc cttacgcacg tctgcggtat ttacaccgc ataatccg
3241 acaccatcga atggtgcaaa accttccgcg gtatggcatg atagcgcgcc gaagagagtc
3301 aattcagggt ggtgaatggt aaaccagtaa cgttatacga tctgcagag tatgccggtg
3361 tctctatca gaccgttcc cgcgtggtga accaggccag ccacgtttct gcgaaaacgc
3421 gggaaaaagt ggaagcggcg atggcgggagc tgaattacat tcccaaccgc gtggcacaac
3481 aactggcggg caaacagtcg ttgctgattg gcgttgccac ctccagtctg gccctgcacg
3541 cgcctgcga aattgtcgcg gcgattaaat ctgcgcccga tcaactgggt gccagcgtgg
3601 tgggtcgtat ggtagaacga agcggcgtcg aagcctgtaa agcggcgggt cacaatctc
3661 tcgcgcaacg cgtcagtggt ctgatcatta actatccgct gtagtaccag gatgccattg
3721 ctgtggaagc tgctgcact aatgtccgg cgttatttct tgatgtctct gaccagacac
3781 ccatcaacag tattatttcc tccatgaag acggtacgcg actggcgtg gagcatctgg
3841 tcgattggg tcaccagcaa atcgcgctgt tagcgggcc attaagtctt gtctcggcgc
3901 gtctcgtct ggctggctgg cataaatatc tcaactgcaa tcaaattcag ccgatagcgg
3961 aacgggaagg cgactggagt gccatgtccg gtttcaaca aaccatgcaa atgctgaatg
4021 agggcatcgt tccactgcg atgctggtg ccaacgatca gatggcgtg ggcgcaatgc
4081 gcgccattac cgatccggg ctgcgcttg gtgcggatat ctcggtagt ggatacgacg
4141 ataccgaaga cagctcatgt tatatcccgc cgtaaccac catcaaacag gattttgcc
4201 tgctggggca aaccagcgtg gaccgcttc tcaactctc tcagggccag gcggtgaagg
4261 gcaatcagct gttgcccgtc tactggtga aaagaaaaac caccctggcg cccaatcgc
4321 aaaccgctc tcccgcgcg ttggccgatt cattaatgca gctggcacga caggttccc
4381 gactggaag cgggcagtga gcgcaacgca ataatgtga gttagtcac tcattaggca
4441 ccccaggctt tacactttat gcttccgct cgtatgtgt gtggaattgt gagcgataa
4501 caattcaca caggaaacag ctatgacct gattacggat tcaactggcg tcttttaca
4561 acgtcgtgac tgggaaaacc ctggcgttac ccaactaat cgccttgca cacaacccc
4621 tttgccagc tggcgaata gcgaagagc cgcaccgat cgccttccc aacagtgcg
4681 cagcctgaat ggcgaatggc gctttgctg gtttccggca ccagaagcgg tcccgaag
4741 ctggctggag tgcgatctc ctgagggcga tactgtcgtc gtcccctcaa actggcagat
4801 gcacggttac gatgcgcca tctacacaa cgtaacctat cccattacgg tcaatccgc
4861 gttgttccc acggagaatc cgacgggtg ttactcgtc acattaatg ttgatgaag
4921 ctggtacag gaaggccaga cgcgaattat tttgatggc gttggaatt

//

Please note: All products are "FOR RESEARCH USE ONLY AND ARE NOT INTENDED FOR DIAGNOSTIC OR THERAPEUTIC USE"

